Bedienungsanleitung Kat. Nr. 38.2047.xx

Bedienungsanleitung Kat. Nr. 38.2047.xx

D

Digitaler Timer und Stoppuhr

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Instruction manual Cat.-No. 38.2047.xx

3. Getting started

- Open the battery compartment using a coin and remove the battery insulation strip.
- 00 MIN and 00 SEC appear on the display.
- Close the battery compartment.
- The device is ready for use.

4. Count Down Timer

- Press the **M** button to adjust the minutes and the **S** button to adjust the seconds (max. 99 min and 59 sec). Press and hold the **M** or **S** button for fast mode.
- Press the **START/STOP** button.
- MIN and SEC will be flashing and the count-down process starts.
- When the timer counted down a beeping alarm will sound (ca. 60 minutes).
- The LED signal lamp flashes.
- Press the **START/STOP** button to stop the alarm.
- The last adjusted alarm time appears on the display.

5. Stopwatch

- Press the **START/STOP** button.
- The count up process starts.

6. Operation

- With the **START/STOP** button you can stop and restart the counting.
- Press the **M** and **S** button at the same time to reset the timer.

7. Attachment

- Use the magnet at the back of the instrument for placement on metal surfaces.

8. Battery replacement

- Change the battery when the display or the alarm tone becomes weaker.
- Make sure the polarities are correct (+pole above).

9. Trouble shooting

Problem

Solution

No display	→ Ensure that the battery polarity is correct → Change the battery
Incorrect indication	→ Change the battery

If your device fails to work despite these measures, contact the retailer where you purchased the product.

11. Specifications

Running time Max. 99 min. 59 sec.

Power consumption Battery 1x CR2032 (included)

Housing dimension 50 x 9 x 80 mm

Weight 29 g (device only)

10. Waste disposal

This product and its packaging have been manufactured using high-grade materials and components which can be recycled and reused. This reduces waste and protects the environment.

Dispose of the packaging in an environmentally friendly manner using the collection systems that have been set up.

Disposal of the electrical device

Remove non-permanently installed batteries and rechargeable batteries from the device and dispose of them separately.

This product is labelled in accordance with the EU Waste Electrical and Electronic Equipment Directive (WEEE).

This product must not be disposed of in ordinary household waste. As a consumer, you are required to take end-of-life devices to a designated collection point for the disposal of electrical and electronic equipment, in order to ensure environmentally compatible disposal. The return service is free of charge. Observe the current regulations in place!

Important information on product safety!

- Do not expose the device to extreme temperatures, vibrations or shocks.
- Protect it from moisture!
- Clean the instrument with a soft damp cloth. Do not use solvents or scouring agents.
- Remove the battery if you do not use the device for a long period of time.

Instruction manual Cat.-No. 38.2047.xx

Disposal of the batteries

Batteries and rechargeable batteries must never be disposed of with household waste. They contain pollutants such as heavy metals, which can be harmful to the environment and human health if disposed of improperly, and valuable raw materials such as iron, zinc, manganese or nickel that can be recovered from waste. As a consumer, you are legally obliged to hand in used batteries and rechargeable batteries for environmentally friendly disposal at retailers or appropriate collection points in accordance with national or local regulations. The return service is free of charge. You can obtain addresses of suitable collection points from your city council or local authority.

The names for the heavy metals contained are:

Cd=cadmium, Hg=mercury, Pb=lead.

Reduce the generation of waste from batteries by using batteries with a longer lifespan or suitable rechargeable batteries. Avoid littering the environment and do not leave batteries or battery-containing electrical and electronic devices lying around carelessly. The separate collection and recycling of batteries make an important contribution to relieving the impact on the environment and avoiding health risks.

5. Stopwatch

- Press the **START/STOP** button.
- The count up process starts.

6. Operation

- With the **START/STOP** button you can stop and restart the counting.
- Press the **M** and **S** button at the same time to reset the timer.

7. Attachment

- Use the magnet at the back of the instrument for placement on metal surfaces.

8. Battery replacement

- Change the battery when the display or the alarm tone becomes weaker.
- Make sure the polarities are correct (+pole above).

9. Troubleshooting

Problem

Solution

No display	→ Ensure that the battery polarity is correct → Change the battery
Incorrect indication	→ Change the battery

If your device fails to work despite these measures, contact the retailer where you purchased the product.

11. Specifications

Running time Max. 99 min. 59 sec.

Power consumption Battery 1x CR2032 (included)

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