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Instruction manual

For your safety

- Keep this instrument and the battery out of reach of children.
- Batteries contain harmful acids and may be hazardous if swallowed. If a battery is swallowed, this can lead to serious internal burns and death within two hours. If you suspect a battery could have been swallowed or otherwise caught in
- the body, seek medical help immediately.

 Batteries must not be thrown into a fire, short-circuited, taken apart or recharged. Risk of explosion!
- Low batteries should be changed as soon as possible to prevent damage caused by leaking. Wear chemical-resistant protective gloves and safety glasses when handling leaking
- Protect it from moisture, extreme temperatures, vibrations

- When the adjusted time is reached, an alarm will ring. The timer is now counting the elapsed time until the alarm is turned off by the START/STOP button. On the display appears automatically the last adjusted time.
- Stopwatch: Press the START/STOP button to start the count up proc
- Press the CLEAR button to clear the display.

Waste disposal

Never dispose of empty batteries rechargeable batteries in household waste, take them to your retail store or to appropriate collection sites.

The symbols for the heavy metals contained are: Cd=cadmium, Hg=mercury, Pb=lead.

Please do not dispose of this instrument in household waste. The user is obligated to take end-of-life devices to a designated collection point for the disposal of electrical and electronic equipment.

Operation

- Open the battery compartment and insert the battery 1,5 V AAA, +/- polarity as indicated. Remove the protective foil from the display.
- Timer: Set the desired time with the button